

Instruction Manual

HI 3826 Ammonia Test Kit for Sea Water



www.hannainst.com

Dear Customer,

Thank you for choosing a Hanna Product.

Please read this instruction manual carefully before using the chemical test kit. It will provide you with the necessary information for correct use of the kit.

Remove the chemical test kit from the packing material and examine it carefully to make sure that no damage has occurred during shipping. If there is any noticeable damage, notify your Dealer or the nearest Hanna office immediately.

Each kit is supplied with:

- 1 beaker (20 mL);
- 1 color comparator cube;
- Reagent 1 (20 mL);
- Reagent 2 (20 mL).

Note: Any damaged or defective item must be returned in its original packing materials.

ISTR3826 12/99 PRINTED IN ITALY

SPECIFICATIONS

Range	0 to 2.5 mg/L NH ₃ -N
Smallest Increment	0.5 mg/L NH ₃ -N
Analysis Method	Colorimetric
Sample Size	10 mL
Number of Tests	25 (average)
Case Dimensions	220x145x55 mm (8.7x5.7x2.1")
Shipping Weight	180 g (6.8 oz.)

SIGNIFICANCE AND USE

The ammonia level in water can vary in nature. Ground water normally contains ammonia due to bacterial decay of plants and animals. However, the presence of ammonia in surface water may be evidence of sanitary pollution due to waste discharges or natural causes.

The Hanna Ammonia Test Kit for Sea Water determines the ammonia concentration in several easy steps. The kit is portable, and can be used in the field as well as in the laboratory.


* mg/L is equivalent to ppm (parts per million)

CHEMICAL REACTION

The ammonia level in mg/L (or ppm), ammonia as nitrogen, is determined by a colorimetric method.

The Nessler reagent reacts with ammonia, under strong alkaline condition, to form a yellow colored complex (see equation below). An addition of a Rochelle salt solution inhibits precipitation of calcium and magnesium ions due to the presence of the alkaline Nessler reagent.

The color intensity of the solution determines the ammonia concentration

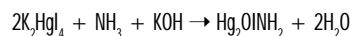


INSTRUCTIONS

READ THE ENTIRE INSTRUCTIONS BEFORE USING THE KIT


- Remove the cap from plastic vessel. Rinse the plastic vessel with water sample and fill to 10 ml mark.




- Add 5 drops of Reagent 2, replace the cap and mix by carefully swirling the vessel in tight circles.

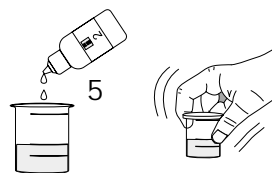


- Add 8 drops of Reagent 1, replace the cap and mix by carefully swirling the vessel.

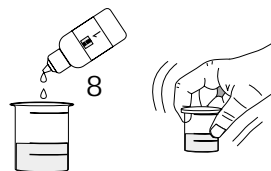

- Remove the cap and transfer the solution into the color comparator cube. Let set for 5 minutes.

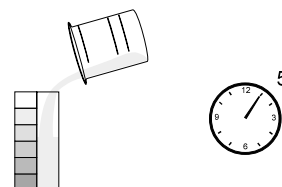

- Determine which color matches the solution in the vessel, and record the results in mg/L (or ppm) NH₃-N

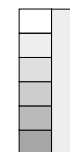

REFERENCES

Standard Methods for the Examination of Water and Wastewater, 16th Edition, 1985, pages 379-382.

HEALTH AND SAFETY

The chemicals contained in this kit may be hazardous if improperly handled. Read Health and Safety Data Sheet before performing this test.