


Instruction Manual

HI 93734

Free & Total Chlorine HR ISM

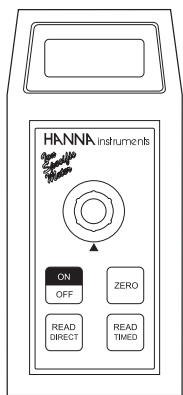

Dear Customer,
Thank you for choosing a Hanna product. This manual will provide you with the necessary information for the correct operation of the meter. Please read it carefully before using the meter. If you need additional technical information, do not hesitate to e-mail us at techserv@hannacan.com. This instrument is in compliance with CE directives EN 50081-1 and EN 50082-1.

PRELIMINARY EXAMINATION

Remove the instrument from the packing material and examine it carefully to make sure that no damage has occurred during shipment. If there is any damage, notify your Dealer.

Each Ion Specific Meter is supplied complete with

- 9V Battery
- Two Sample Cuvets and Caps
- One Transport Cap

Note: Conserve all packing material until the instrument has been observed to function correctly. Any defective item must be returned in its original packing.

GENERAL DESCRIPTION

The HI 93734 meter measures the Free and Total Chlorine (Cl_2) content in water and wastewater.


The meter uses an exclusive positive-locking system to ensure that the cuvet is in the same place every time it is placed into the measurement cell.

The reagents are in liquid form and are supplied in bottles. The amount of reagent is precisely dosed to ensure maximum repeatability.

Display codes aid the user in routine operations.

The meters have an auto-shut off feature that will turn the instrument off after 10 minutes of non-use.

SPECIFICATIONS

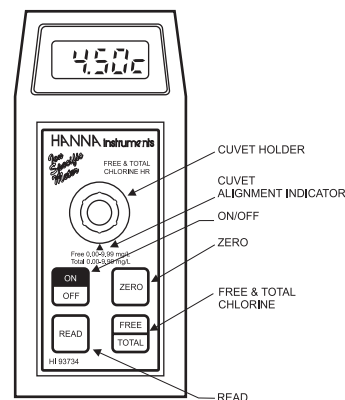

SPECIFICATIONS

Range	Free Cl_2 0.00 to 9.99 mg/L Total Cl_2 0.00 to 9.99 mg/L
Resolution	0.01 mg/L
Accuracy	± 0.03 mg/L $\pm 3\%$ of reading
Typical EMC	± 0.01 mg/L
Deviation	
Light Source	Light Emitting Diode @ 555 nm
Method	Adaptation of the EPA recommended DPD method 330.5. The reaction between the chlorine and the DPD reagent causes a pink tint in the sample
Light Detector	Silicon Photocell
Environment	0 to 50°C (32 to 122°F); max 95% RH non-condensing
Battery Type/Life	1 x 9 volt/40 hours
Auto-Shut off	After 10' of non-use
Dimensions	180 x 83 x 46 mm (7.1 x 3.3 x 1.8")
Weight	290 g (10 oz.).

REQUIRED REAGENTS

Code	Description	Quantity
HI 93734A-0	Free & Total Chlorine DPD 3 drops Reagent	
HI 93734B-0	Free & Total Chlorine Reagent	5 mL
HI 93734C-0	Total Chlorine only	1 drop

REAGENT SETS

HI 93734-01	Reagents for 100 Free & Total Chlorine tests
HI 93734-03	Reagents for 300 Free & Total Chlorine tests

DISPLAY CODE GUIDE

---	This indicates that the meter is in a ready state and zeroing can be performed.
5 1 P	Sampling in Progress. This prompt appears each time the meter is performing a measurement.
-00-	This indicates that the meter is in a zeroed state and measurement can be performed. A zero reading was not taken. Insert a sample before adding reagent and press ZERO.
2 ER 0	Under range. A blinking "0.00" indicates that the sample absorbs less light than the zero reference. Check the procedure and make sure you use the same cuvet for reference (zero) and measurement.
0.00	Over range. A flashing value higher than the maximum concentration readable (see specifications) indicates that the sample absorbs too much light, meaning that the concentration is too high. Dilute the sample.
330	Light over range. The cuvet is not inserted correctly and an excess ambient light is reaching the detector. If the cover is properly installed, then contact your dealer or the nearest Hanna Customer Service Center.
CAP	Light under range. The zero sample is too dark for proper zeroing. If this is not the case, contact your dealer or the nearest Hanna Customer Service Center.
L 0	The "V" indicates that the battery voltage is getting low and the battery needs to be replaced.
V 250	This indicates that the battery is dead and must be replaced.
-88-	Note: once this indication is displayed, the meter will lockup. Change the battery to restart.

HANNA
instruments

www.hannacan.com



This Instrument is in Compliance with the CE Directives

WARRANTY

HI 93734 is warranted for two years against defects in workmanship and materials when used for its intended purpose and maintained according to instructions.

This warranty is limited to repair or replacement free of charge.

Damages due to accident, misuse, tampering or lack of prescribed maintenance are not covered.

If service is required, contact the dealer from whom you purchased the instrument. If under warranty, report the model number, date of purchase, serial number and the nature of the failure. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the instrument is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization Number from the Customer Service department and then send it with shipment costs prepaid. When shipping any instrument, make sure it is properly packaged for complete protection.

To validate your warranty, fill out and return the enclosed warranty card within 14 days from the date of purchase.

All rights are reserved. Reproduction in whole or in part is prohibited without the written consent of the copyright owner, Hanna Instruments Inc.

Hanna Instruments reserves the right to modify the design, construction and appearance of its products without advance notice.

OPERATIONAL GUIDE

MEASUREMENT PROCEDURE

- Turn the meter on by pressing ON/OFF.



- When the LCD displays "-- --", it is ready.




- Select Free or Total Chlorine by pressing FREE/TOTAL. An "I" or a "C" will appear on the right corner to indicate Free or total chlorine, respectively.



- Fill the cuvet with 10 mL of unreacted sample, up to the mark, and replace the cap.




- Place the cuvet into the holder and ensure that the notch on the cap is positioned securely into the groove.

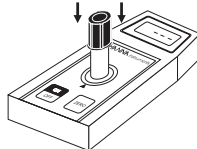

- Press ZERO and "SIP" will appear on the display.




- Wait for a few seconds and the display will show "-0.0-". Now the meter is zeroed and ready for measurement.

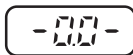

- Remove the cuvet.



- To measure Free Chlorine, add to the other cuvet 3 drops of HI 93734A-0 and 5 mL of HI 93734B-0, using the 5 mL syringe. To measure Total Chlorine,

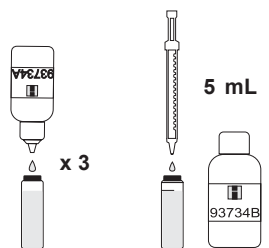

add also 1 drop of HI 93734C-0. Shake for 10 seconds.



Note: To measure exactly 5 mL of reagent with the syringe, push the plunger completely into the syringe and insert the tip into HI 93734B-0 reagent bottle. Pull the plunger out until the lower edge of the seal is on the 5 mL mark of the syringe.

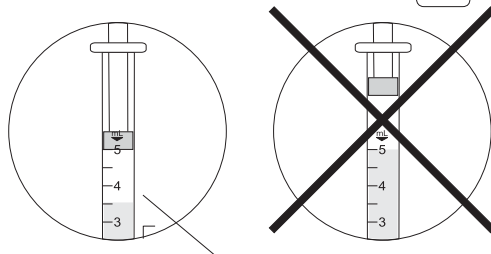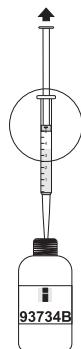


probable level of liquid taken up by syringe

- Add 5 mL of unreacted sample, up to the 10 mL mark of the cuvet, by means of the plastic pipette. Replace the cap and shake gently again.

5 mL of sample



- Reinsert the cuvet into the instrument.

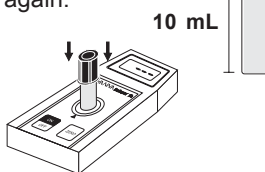

- Wait for approximately 30 seconds and then press READ. The display will show "SIP" during measurement.



- The instrument directly displays concentration in mg/L of Free or Total chlorine on the Liquid Crystal Display.

Notes: Free and total chlorine have to be

measured separately with fresh unreacted samples following the above procedure if both values are requested.

Rinse the plastic pipette 2 or 3 times with sample before adding it to the cuvet.

INTERFERENCES

Interference may be caused by:

Bromine
Iodine
Fluorine
Ozone

Oxidized manganese and Chromium Alkalinity above 250 mg/L or acidity above 150 mg/L will not reliably develop the full amount of color or it may rapidly fade. To resolve this, neutralize the sample with diluted HCl or NaOH.

In case of water with hardness greater than 500 mg/L CaCO₃, shake the sample for approximately 1 minute after adding the reagents.

ACCESSORIES

- 5 mL syringe
- 1 plastic pipette

TIPS FOR AN ACCURATE MEASUREMENT

The instruction listed below should be carefully followed during testing to ensure best accuracy.

- Do not touch the cuvet walls with hands.
- In order to maintain the same conditions during the zeroing and the measuring phases, it is necessary to close the cuvet to prevent any contamination.
- Do not let the test sample stand too long after reagent is added or accuracy will be lost.
- Whenever the cuvet is placed into the measurement cell, it must be completely free of fingerprints, oil or dirt. Wipe it thoroughly with HI 731318 or a lint-free cloth prior to insertion.
- It is important that the sample does not contain any debris. This would corrupt the readings.
- It is possible to take multiple readings

in a row, but it is recommended that a zero reading be taken for each sample and that the same cuvet is used for zeroing and measurement.

- It is important to discard the sample immediately after the reading is taken because the glass might become permanently stained.
- Shaking the cuvet can generate bubbles in the sample, causing higher readings. To obtain accurate measurements, remove such bubbles by swirling or by gently tapping the vial.
- All the reaction times reported in this manual are referred to 20°C (68°F). As a general rule of thumb, they should be doubled at 10°C (50°F) and halved at 30°C (86°F).

ACCESSORIES

REAGENT SETS

HI 93734-01 Reagents for 100 Free & Total Chlorine tests
HI 93734-03 Reagents for 300 Free & Total Chlorine tests

OTHER ACCESSORIES

HI 710009 Blue rubber boot
HI 710010 Orange rubber boot
HI 721310 9V battery (10 pcs)
HI 731318 Tissue for wiping cuvetts (4 pcs)
HI 731321 Glass cuvetts (4 pcs)
HI 731325 Caps for cuvetts (4 pcs)
HI 93703-50 Cuvetts cleaning solution (230 mL).

CE DECLARATION OF CONFORMITY

Recommendations for Users
Before using these products, make sure that they are entirely suitable for the environment in which they are used. Operation of these instruments in residential area could cause unacceptable interferences to radio and TV equipments, requiring the operator to take all necessary steps to correct interferences. Any variation introduced by the user to the supplied equipment may degrade the instruments' EMC performance. To avoid damages or burns, do not perform any measurement in microwave ovens.

